QsepSeries *Fragment Analyzer*

DNA/RNA Samples

Gel Preparation

Manual Loading

Electrophoresis

Gel Imaging

All in One Step

Qsep Series DNA/RNA Fragment Analyzer

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Routine Applications

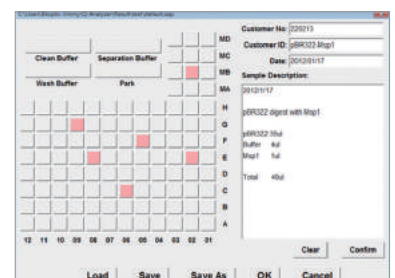
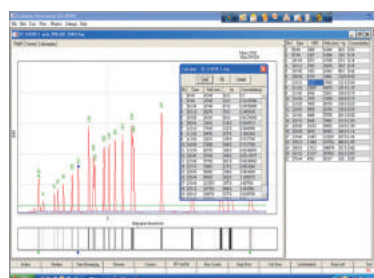
- PCR Product Screening
- Genotyping
- Next Generation Sequencing (NGS) QC
- Restriction Fragment Length Polymorphism (RFLP)
- Oligonucleotides Analysis
- Genomic DNA Analysis
- Plasmid Purification & Vector Cloning Analysis
- RNA Analysis
- Protein Profiling (fluorescence detection)

Specific Applications

- CRISPR technology
- cfDNA (Circulating Cell Free DNA)
- Chromatin immunoprecipitation (ChIP)
- Glycan analysis (N-Glycan)
- Immuno Assay : Ag and Ab Reaction
- Genetic profiling /disease screening
- Pathogen detection/typing (HBV, HCV, HPV, TB, AIV, STD etc.)
- Agriculture product detection (GMO, 한우판정, 식물품종 판별)
- Food Safety analysis
- SNP screening
- Microsatellite test (STR, SSR, MSI)
- 3-bp repeat disease (Fragile X, HD, SCA)

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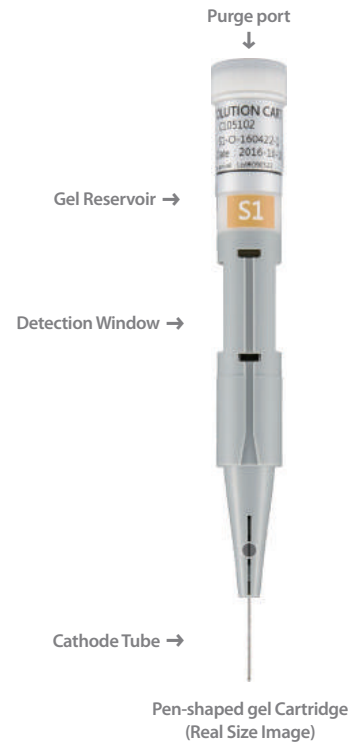
4-Channel (Qsep400)

Cartridge Selection Guide

Cartridge	Standard (S2)	High Resolution (S1)	Kilo Base (S3)	Fast (F3)	High Sensitivity (N1)	RNA (R1)	Protein (P2)
Sample size range	10-15000 bp	10-15000 bp	10-23000 bp	10-5000 bp	10-5000 bp	N/A	14-150 kDa
L.O.D.	0.1 ng/μl*	0.1 ng/μl*	0.1 ng/μl*	0.1 ng/μl*	5 pg/μl*	5 ng/μl	0.5 ng/ul
Resolution	4-10 bp	1-4 bp	10-50 bp	≥ 50 bp	≥ 10 bp	N/A	N/A
Analysis Time (per sample)	2-3min	3-5min	5-8min	1-2min	3-5min	5-10min	6-10min

L.O.D.: 2pg/μl (If diluted in the water) * Determined by the DNA ladder (15-622 DNA size marker : C109200) as sample
 KIT 구성품 : Alignment marker, Separation buffer, Dilution buffer, Mineral oil, buffer tray, PCR vial and dropper

Application	Standard (S2)	High Resolution (S1)	Kilo Base (S3)	Fast (F3)	High Sensitivity (N1)	RNA (R1)	Protein (P2)
PCR product screening	●	●		●			
RFLP	●	●		●			
SNP		●					
Plasmid purification & Vector cloning analysis			●				
Library & Fragmentation	●	#					
Genomic DNA analysis	@		●				
Cell free DNA detection					●		
RNA Analysis						●	
Low concentration sample					●		
CRISPR QC & Analysis	●	●				●	
Large size fragment analysis			●				
Replace SDS-PAGE							●
IgG Purity Test							●

#. Higher accuracy, but longer separation time. @. Genomic DNA size below 15kb

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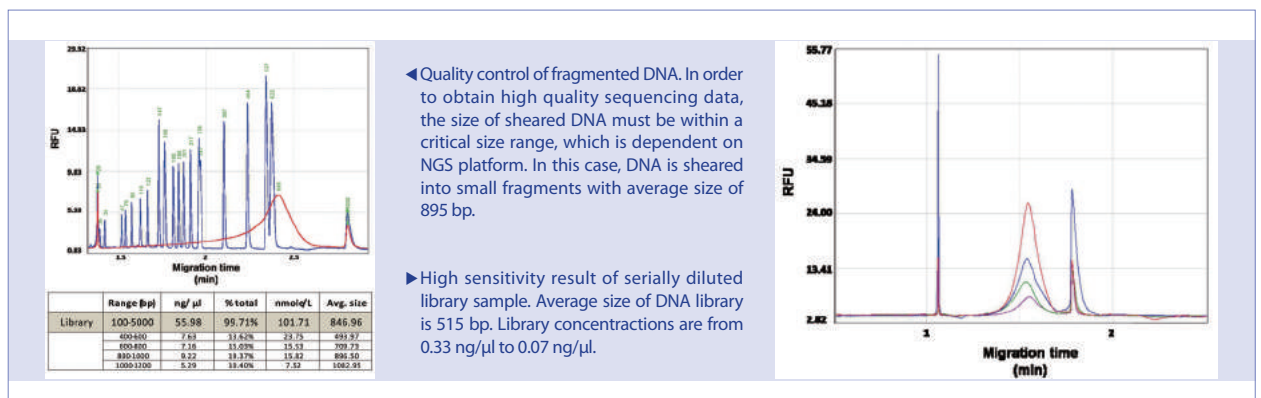
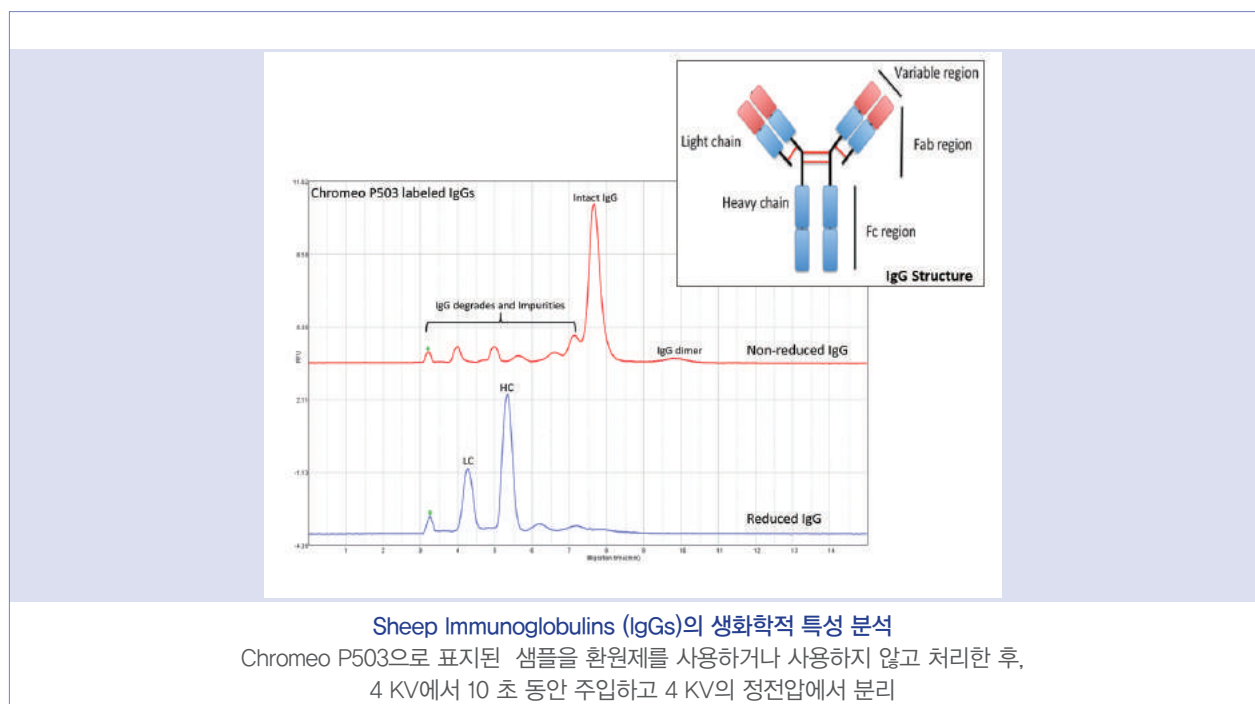
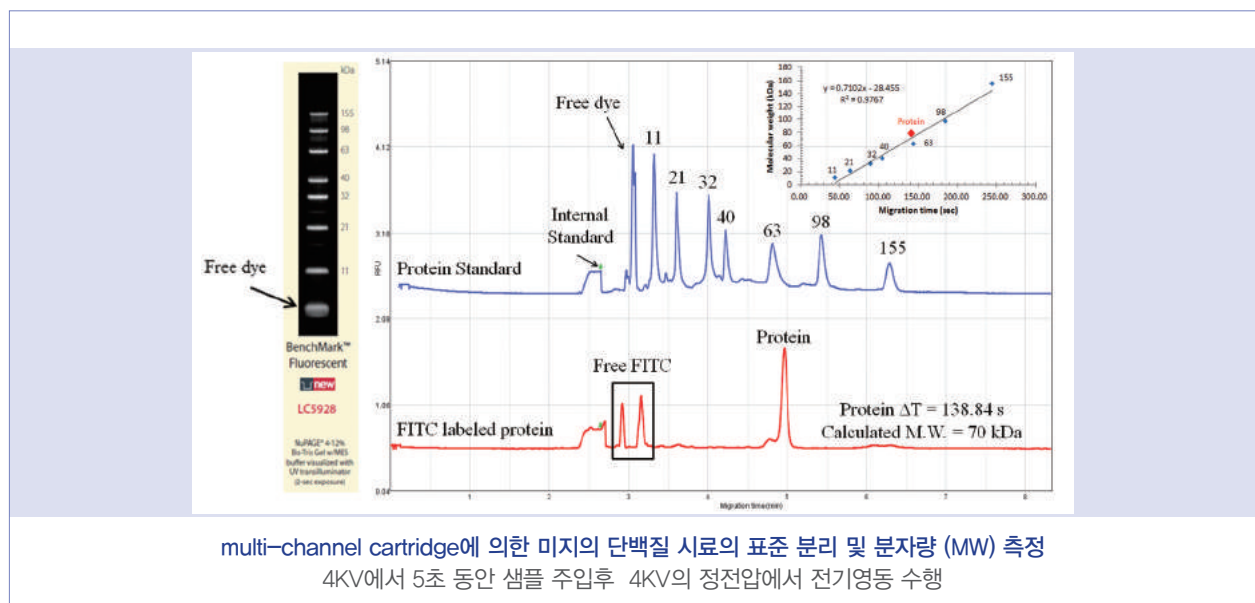
◀Representative cell-free DNA sample result using the Qsep100 analyzer with N1 high sensitivity cartridge kit. Low amounts cfDNA, yields less than 5 pg/μl, is directly detectable by Qsep100 instrument.

Fluorescence Detection of Protein

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Bi_optic

Highlights

	Qsep1	Qsep100	Qsep400
System Type	1 Channel Portable System	1 Channel Standard System	4-Channel System
Automated Sampling	1~8 samples	1~96 samples	1~96 samples
Disposable Gel-Cartridge*	Single Channel: S1, S2, N1,R1	Single Channel: S1, S2, F3, N1,R1	Four Channel: S1, S2, F3, N1,R1
Rapid Analysis	2~7 min/sample	2~7 min/sample	2~7 min/4 samples
Resolution	1~4bp (Between 100~500bp)	1~4bp (Between 100~500bp)	1~4bp (Between 100~500bp)
Sensitivity	5pg/μl	5pg/μl	5pg/μl
Minimum Sample Volume	1 μl (Micro-Vial:C104250) 20 μl (standard PCR tube)	1 μl (Micro-Vial:C104250) 20 μl (standard PCR tube)	1 μl (Micro-Vial:C104250) 20 μl (standard PCR tube)
Sample Consumption	≤1pl	≤1pl	≤1pl

* (100 to 300 sample capacity without gel preparation)

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